# SOEHNLE

## PROFESSIONAL

|                                |    |
|--------------------------------|----|
| Bedienungsanleitung .....      | 2  |
| Operating Instructions .....   | 5  |
| Mode d'emploi .....            | 8  |
| Gebruiksaanwijzing .....       | 11 |
| Instrucciones de manejo .....  | 14 |
| Istruzioni d'uso .....         | 17 |
| Návod k obsluze .....          | 20 |
| Käyttöohje .....               | 23 |
| Bruksanvisning (S) .....       | 26 |
| Bruksanvisning (N) .....       | 29 |
| Brugsanvisning .....           | 32 |
| Οδηγίες χρήσης .....           | 35 |
| Instrukcja obsługi .....       | 38 |
| Kullanma Kılavuzu .....        | 41 |
| Manual de<br>instruções .....  | 44 |
| Használati<br>utasítás .....   | 47 |
| Lietotāšanas instrukcija ..... | 50 |
| Naudojimo instrukcija .....    | 53 |
| Kasutusjuhend .....            | 56 |
| Návod na obsluhu .....         | 59 |
| Navodila za uporabo .....      | 62 |

7755 • 7755 plus • 7756 • 7757

## Montageanleitung 7755 plus

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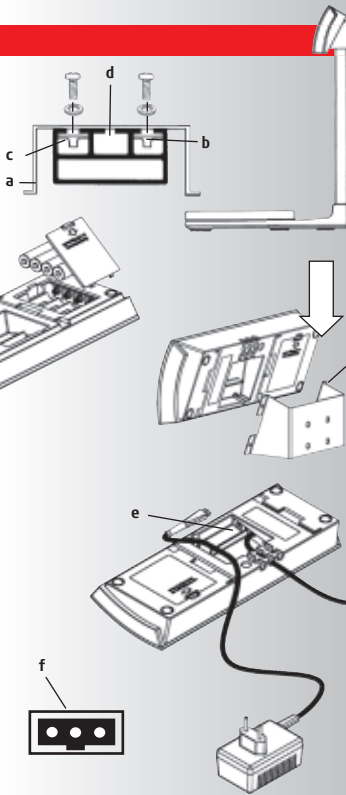
## Installation instructions for 7755 plus

Insert the screws into the holes in the display holder (a) and hand-tighten them using the enclosed, self-locking lock nuts (b). Run the measuring cable up the rear side of the display stand (d). Insert the display holder (a) with the locking nuts (b) into the guide groove (c) of the display stand and fasten it. Push the display unit down into the display holder (a) until it locks into place.

### Getting started

Insert batteries (4 x 1.5 V AA) or connect the unit to a power supply. Position the scale at its place of use. Make sure it is horizontal. The scale is ready for operation.

If you want to fix the display unit to a wall, attach the enclosed display holder (a) to the wall and push the display unit down into it until it locks into place.

### Plug-in power supply unit

The socket (f) for connecting the supplied plug-in power supply unit is located at the base of the display unit (e). Insert the cable and run it into the cable grip. Always use an original Soehnle Professional power supply unit. The power supply unit is not waterproof and therefore must not come into contact with water.

Should the power supply unit become wet or show visible signs of damage, do not operate the scale with the power supply unit. Please contact your dealer or service partner.

### Power supply

You can also operate the scale using Soehnle Professional rechargeable batteries besides the plug-in power supply unit and zinc/carbon or alkali/manganese batteries (4 x 1.5 V AA).

Please note that the batteries are not charged when the power supply unit is switched on. You require a separate battery recharging unit.

### Recharging kit (optional)

You can obtain Soehnle Professional rechargeable batteries as part of a separate kit, consisting of a recharging unit, Soehnle Professional rechargeable batteries and a safety sticker. When inserted, Soehnle Professional rechargeable batteries have sufficient charge for up to 200 hours of weighing (depending on the recharging level of the batteries). During recharging (approx. 15 hrs), an indicator lamp is lit on the recharging unit. Start a new loading cycle by removing and then inserting the flat plug on the display unit. If you do not use the rechargeable batteries for a long time, they will lose some of their capacity. The batteries are fully charged after approx. 3 to 5 recharging cycles.

Replacement parts: Soehnle Professional rechargeable batteries, recharging unit

When you insert rechargeable batteries into the unit, attach the supplied sticker to the battery compartment.

1. The sticker reads as follows: Caution! Rechargeable Soehnle Professional batteries in use, do not replace with other batteries.

**Caution:** if you use normal batteries with the battery recharging unit, there is a danger of leakage and explosion. If you start using normal batteries again, replace the battery recharging unit with the plug-in power supply unit supplied with the scale.

2. The power supply unit is a battery recharging unit and can be recognised by the LED on the housing.

#### Technical data

Refer to the information on the display unit for weight range and numeric intervals.

The power supply data is given on the plug-in power supply unit.

Working temperature: 5 - 35°C

Storage temperature: -10 - 55°C

Dispose of used (rechargeable) batteries in accordance with the national and local regulations for environmental protection and recycling.

**Weighing**

Switch the scale on.  
The test display "888888" appears. Then "o" appears.  
Scale is ready for weighing.

**Tare weighing and add-on weighing**

Place the container on the base. Press the  key. "o" is displayed. Place the weighing goods on the scale.  
Add-on weighing: Press the  key again until "o" is displayed. Add the additional weighing goods. You can keep doing this up to the maximum load weight. After the first tare weighing, an "N" (net) appears in the display for information purposes. "N" is only cleared when you switch the scale off.

**Check weighing**

Place the reference object on the scale. Tare. Remove the reference object. Place the object under test on the scale and read off the plus/minus difference on the display.

**Total weight display**

When objects are on the scale, press the  key. "o" is displayed. Remove the goods for weighing. The total weight appears with a minus sign.

**Switching off**

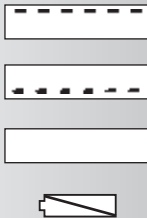
When the display shows "o", press the  key. When operating the scale with batteries or rechargeable batteries, the auto-off function is active. When the display shows "o", the scale switches off automatically after 2 minutes.

**Changing units**


This is possible at any time. Press the  key to change units. The last unit used remains active when you switch the scale off and appears again the next time you switch the unit on.

**Fault messages**


- **Overload:** please observe the weight range of the scale (see name-plate).
- **Underload:** switch the scale on again. If this does not solve the problem, contact your dealer or service partner.
- **With an unclear display** (see example on left): switch off the power supply to the scale for 10 seconds by removing the (rechargeable) batteries and/or pulling out the power plug.
- **(Rechargeable) batteries empty:** Please replace or charge them (4 x 1.5 V AA). Please make sure that you do not try to charge the rechargeable batteries with the power supply unit plugged in. You require a separate recharging unit. This is available as an accessory.

**Important additional information accu set 2557**

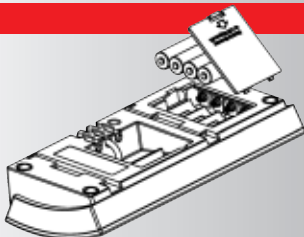**Attention:** The accu set is only pre-loaded. The scale can be put into operation immediately. Before the first usage we advice to load the accu for at least 4 hours in order to achieve the full accu capacity.  
**If the appliance is operated with the charge adapter, only use rechargeable batteries. Normal batteries will destroy the appliance.**

## User calibration

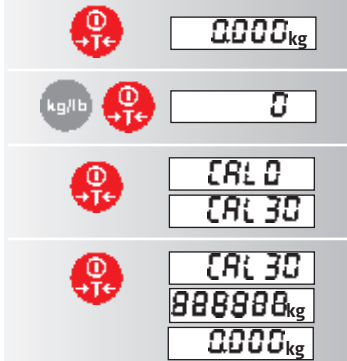If necessary, you can readjust the scale. To do this, you require a calibration weight of a weight equal to the maximum weight range.

1. Switch the scale on until "o" is displayed.
2. Press both keys (0 + T) for 20 seconds. There must be NO weight on the scale. "o" is displayed.
3. Press the 0 key. "CAL o" appears and then the maximum weight range of the scale.
4. Place the calibration weight on the scale. This weight must be equal to the maximum weight range. Press the 0 key. "CAL XX" appears (XX = maximum weight range), then "888888" and finally "o". Then, take the weight from the scale and switch the scale off.

User calibration is always displayed in kg. Press the T key to cancel the operation.

## Cleaning

It is sufficient to clean the scale using a damp cloth and a normal cleaning agent. Do not use any scourers.

## Maintenance

The scale does not require any routine maintenance. However, we recommend to check the scale's accuracy at regular intervals. The regularity of these checks is dependent on the level of use and the state of the scale. If any inaccuracies occur, please contact your dealer or the Soehnle Professional Service Department.

## Note

The display value may be influenced by extreme electrostatic and electro-magnetic interference, e.g. operation of a radio unit or a mobile phone in close vicinity to the unit. When the electrostatic interference is no longer present, the unit can again be used for its intended purpose. You may need to switch it on again.

In addition, air currents, vibrations, rapid changes in temperature and direct sunlight may have an effect on the accuracy of the scale. The scale is protected to IP 54 protection class. Therefore, avoid high levels of humidity, vapours, liquids and strong contamination. These influences may damage the measuring components and render your warranty claim null and void.

**Interference suppression of the unit corresponds to the valid EU directive 89/336/EEG.**

## Warranty - Liability

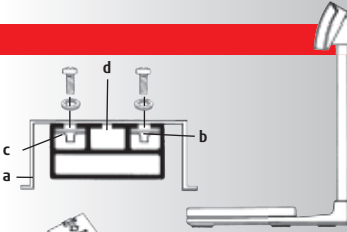
If a fault or defect is present on receipt of the unit which is within Soehnle Professional's scope of responsibility, Soehnle Professional shall be obliged to repair the fault or supply a replacement unit. Replaced parts shall be the property of Soehnle Professional. Should the fault repairs or replacement delivery not be successful, the statutory provisions shall be valid. The period of warranty shall be 2 years, beginning on the day of purchase. Please retain your receipt as a proof of purchase. Should your scale require servicing, please contact your dealer or Soehnle Professional Customer Services.

### No responsibility shall be accepted for damage caused through any of the following reasons:

Unsuitable or improper storage or use, incorrect installation or commissioning by the owner or third parties, natural wear, changes or modifications, incorrect or negligent handling, overuse, chemical, electrochemical or electrical interference or humidity, unless they can be traced back to Soehnle Professional.

If operating, climatic or any other influences lead to a major change in conditions or material quality, the warranty for perfect unit functioning shall be rendered null and void. If Soehnle Professional provides an individual warranty, this means that the unit supplied will be free of faults for the length of the warranty period.

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**Soehnle Professional GmbH & Co. KG**